
MOLD RELEASE

PURPOSE: A final rinse for the metal embedding molds. After staining, paraffin will not stick to the molds.

PRINCIPLE: The glycerin in the mold release coats the metal molds.

SOLUTION:

Mold Release

70-95% Alcohol 4000.0 ml
Glycerin 50.0 ml
Mix well, cover. Stable for 2
months.

CAUTION: Flammable.

SAFETY /PPE: Wear gloves, goggles, and lab coat. Avoid contact and inhalation.

Prepared: _____ By: _____

Approved: _____ By: _____

MOLD RELEASE

70-95% Alcohol 4000.0 ml
Glycerin 50.0 ml

Mix well, cover. Stable for 2 months.

CAUTION: Flammable.

DATE: _____

TECH: _____

EXPIRATION: _____